

For more information



Rebuilding the Cell-Free System and the Applications for R&D of Biologics


Reconstituted cell-free protein synthesis kit

PUREfrex[®]

in vitro protein selection technology

PUREfrex[®]RD

Takashi Ebihara, Ph.D.
COO
GeneFrontier Corporation

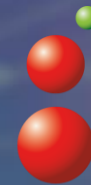


GeneFrontier

PEGS Europe
14-16 Nov 2022

PUREfres[®]

-Customize expression toolbox for your research-

Only necessary molecules
for transcription/translation

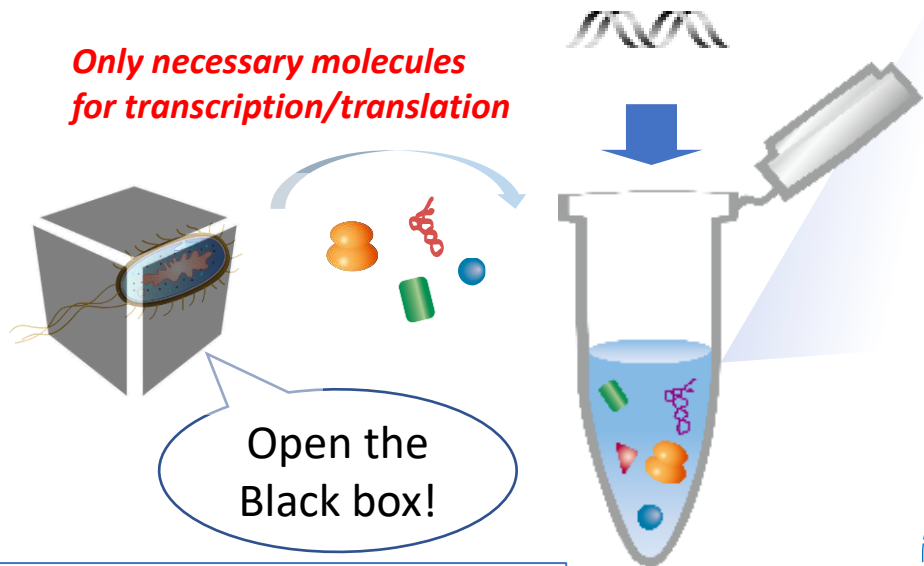


Open the
Black box!




- ✓ Rebuilt cell-free system
- ✓ Tunable for your biologics
- ✓ Simple, Fast
- ✓ Suited for High throughput system
- ✓ Very low contaminants



Totally constructive, molecular based system

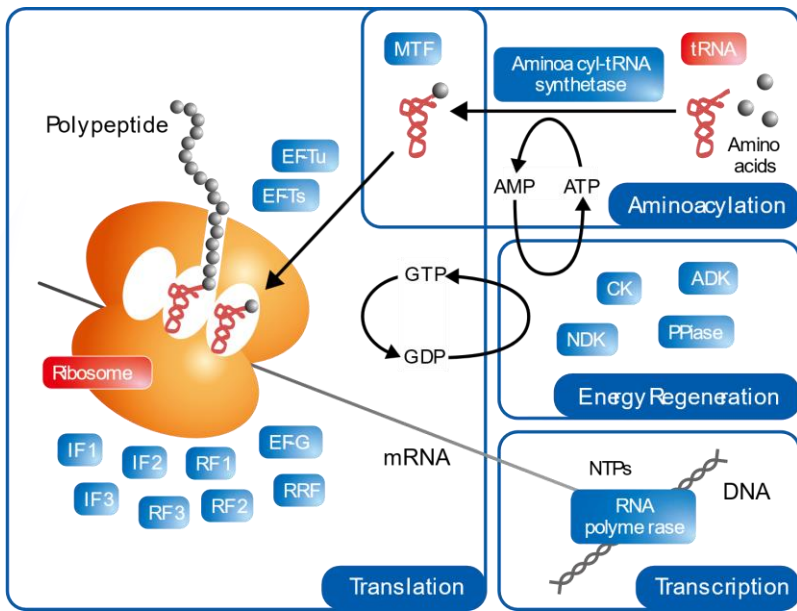

For more information

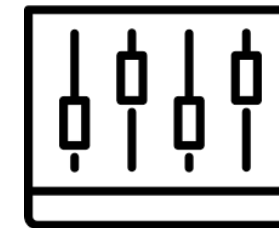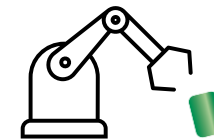
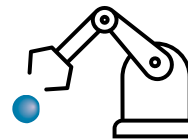


Translation factors
Chaperones
Detergent
Temperature/pH

<Ex, Non-natural AA introduction>

Translation - RF1


Suppressor tRNA + non-natural amino acid

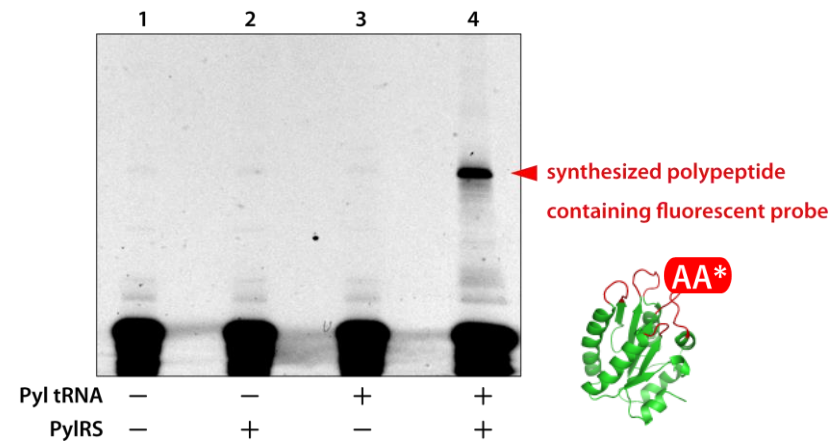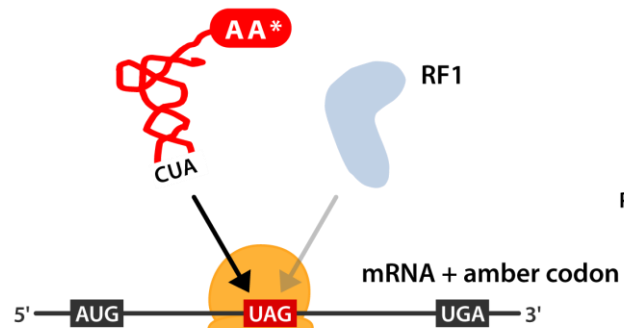

PUREfres[®]

-Customize expression toolbox for your research-


Only necessary molecules for transcription/translation

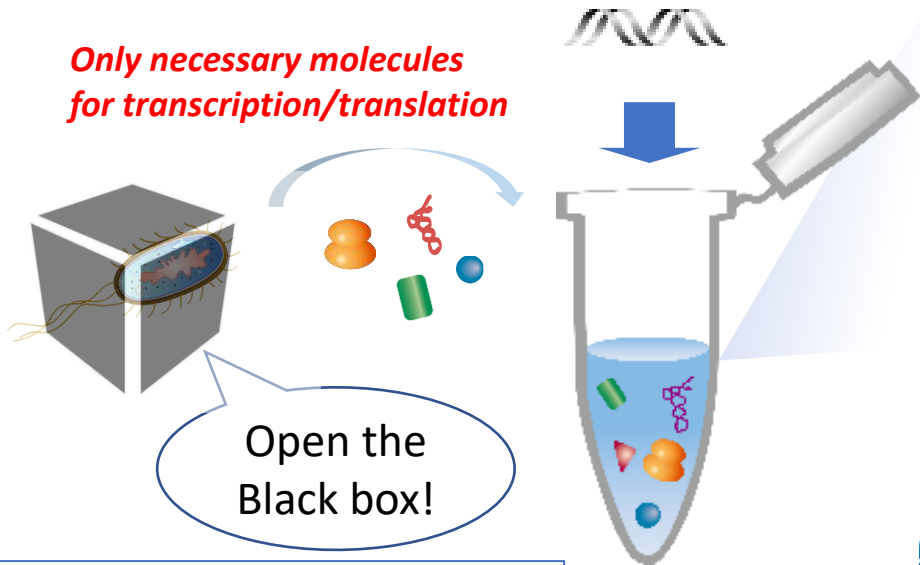




Open the Black box!

- ✓ Rebuilt cell-free system
- ✓ Tunable for your biologics
- ✓ Simple, Fast
- ✓ Suited for High throughput system
- ✓ Very low contaminants

Totally constructive, molecular based system

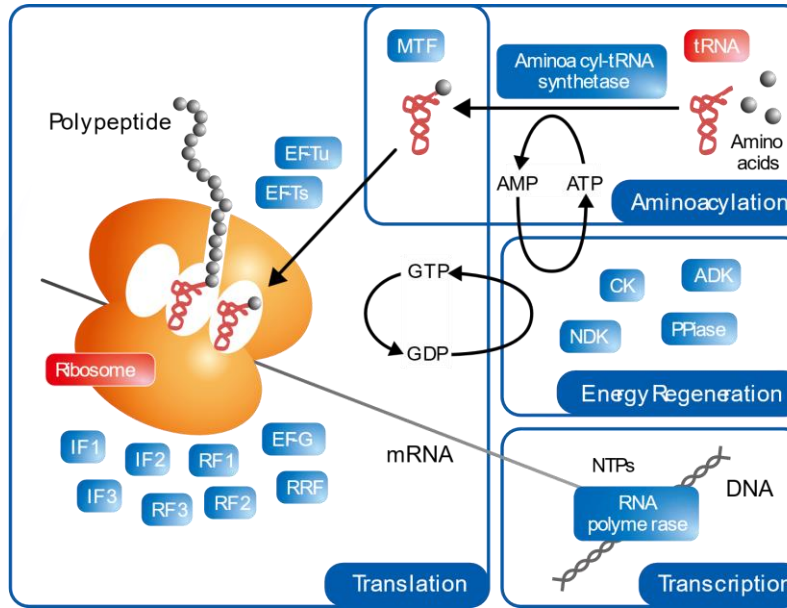


For more information

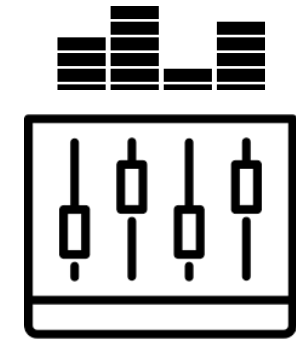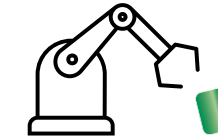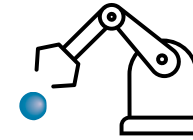

Translation factors
Chaperones
Detergent
Temperature/pH


<Ex, Membrane protein with Nanodisc; artificial membrane-like structure>

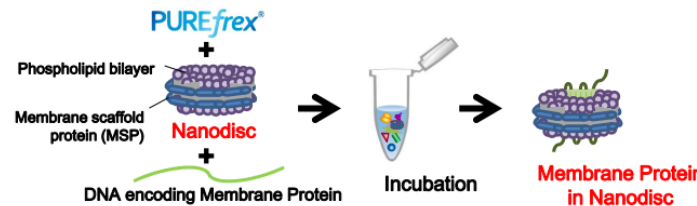

The condition of membrane protein synthesis

Reaction mix	Template DNA	Incubation
PUREfres [®] 2.0 +Nanodisc (MSP1E3D1-His POPC*, final 10 μM)	PCR product	37°C, 4 h

*Ref: Denisov et al. (2007) *J.Biol.Chem.*, vol. 282, p. 7066.

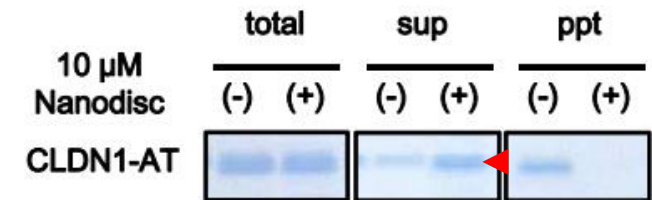

Solubilized hCLDN1 was synthesized using PUREfres[®] and Nanodisc.


Poster MBSJ 2016



PUREfres[®]

-Customize expression toolbox for your research-

Only necessary molecules for transcription/translation


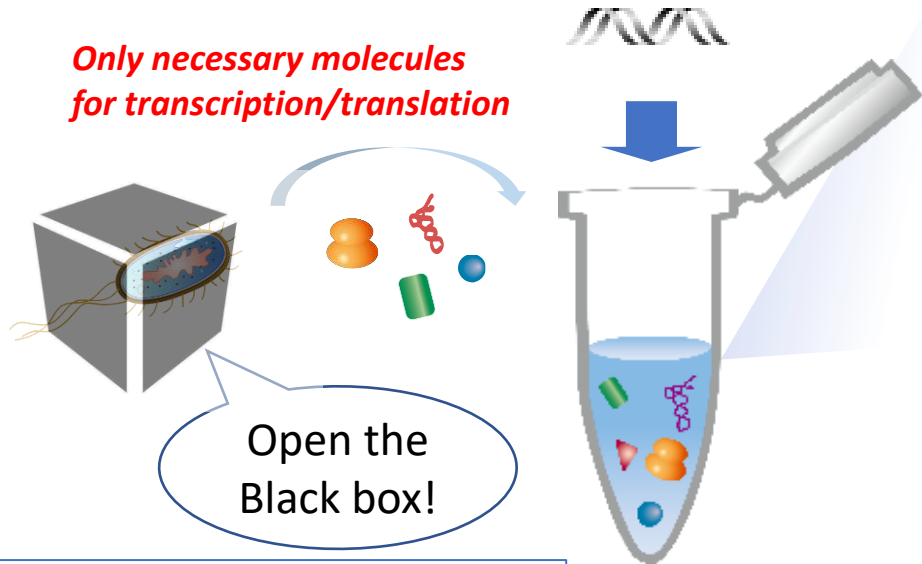


Open the Black box!

- ✓ Rebuilt cell-free system
- ✓ Tunable for your biologics
- ✓ Simple, Fast
- ✓ Suited for High throughput system
- ✓ Very low contaminants



Totally constructive, molecular based system

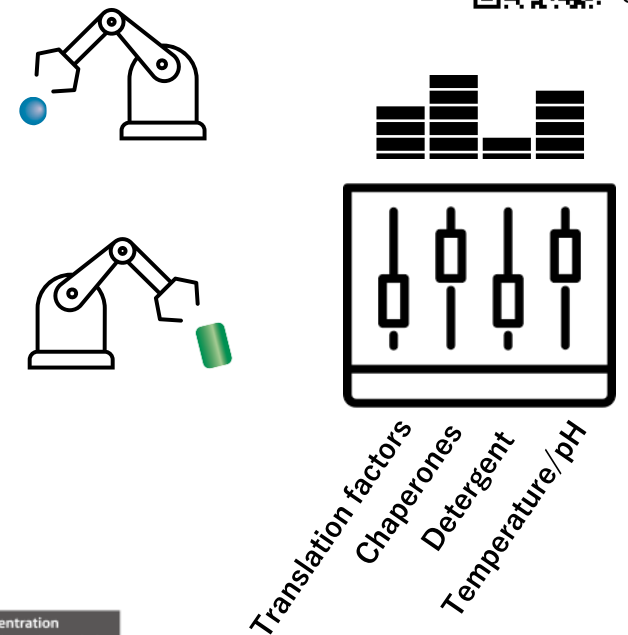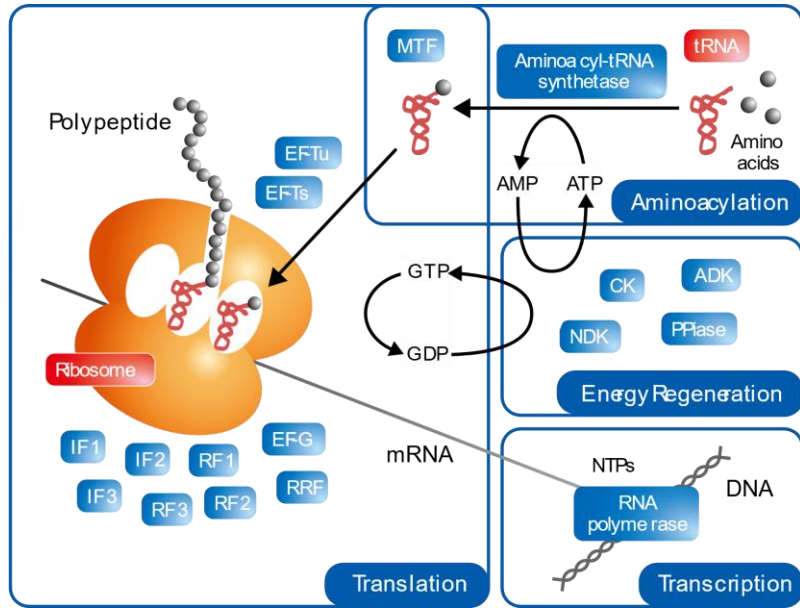

For more information 




Experimental conditions for protein synthesis

Reaction mixture	Incubation	Template DNA
PUREfres [®] 2.1 (4 mM GSH) + Sufractants	37°C 4 h	sfGFP PCR product (1 ng/μL)


→ Measurement of GFP fluorescence

- Most surfactants did not inhibit the protein synthesis reaction by PUREfres[®] below the CMC.
- Some surfactants such as Triton X-100 and Tween 20 could be used even above the CMC.



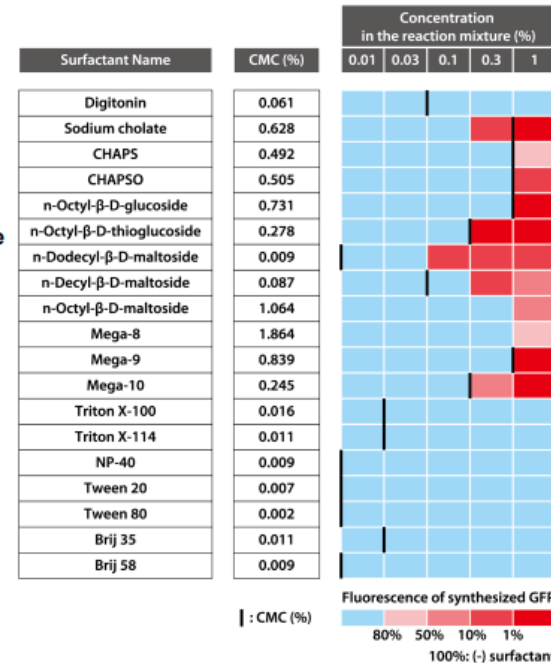




-Having good productivity-

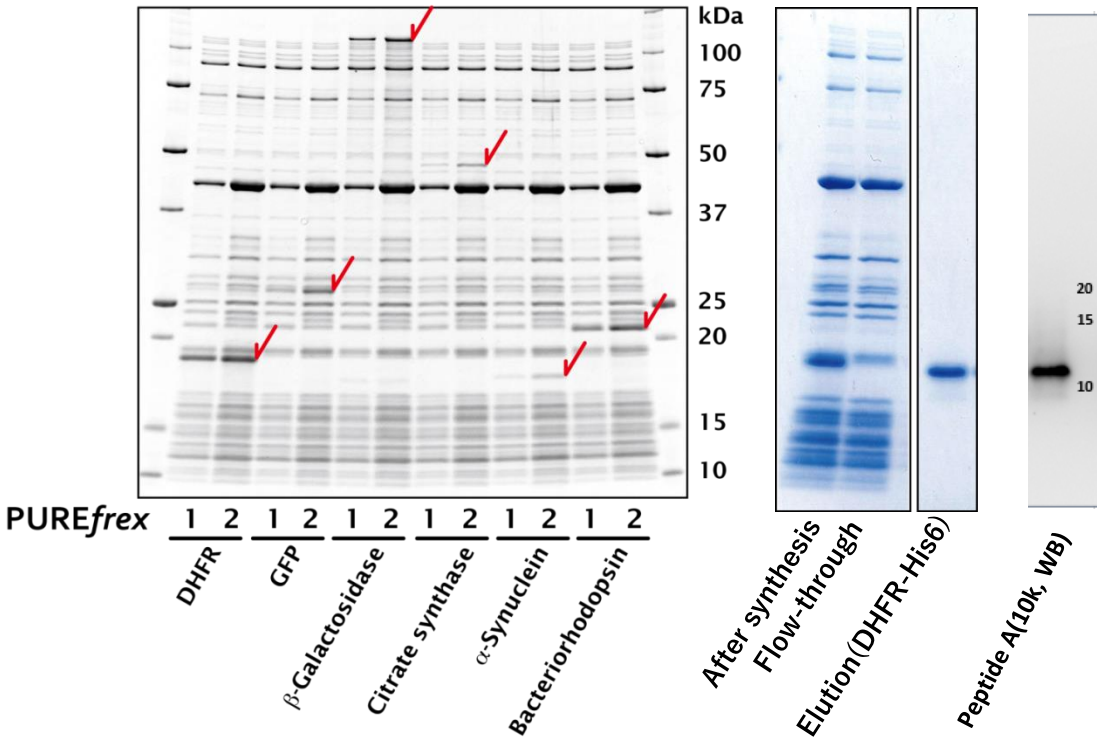

- Reaction at 37°C for 4 h
- 0.5 μ L of reaction mix/ lane
- stained with Oriole (Bio-Rad) and analyzed with an image analyzer (LAS)

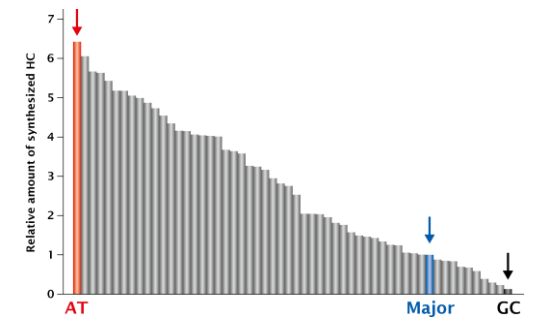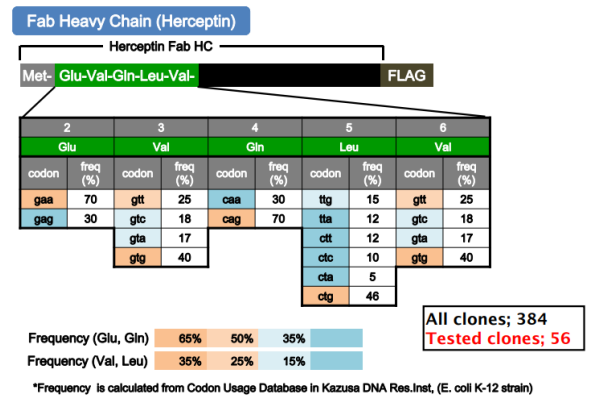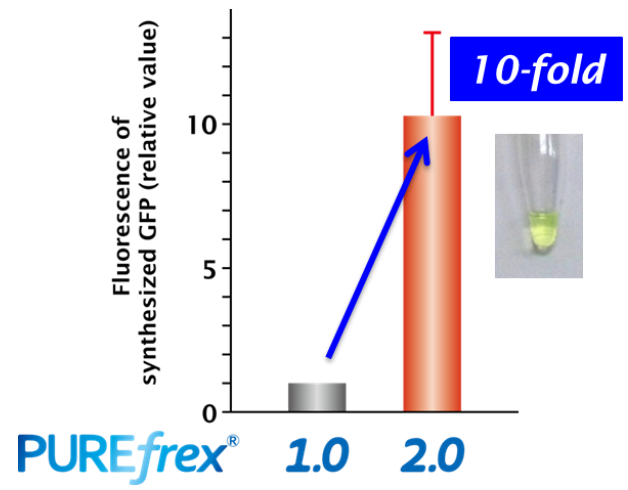

Design of DNA template is important.
Manual is Free to download from our Web site.





- ✓ Good expression for many proteins, small to large.
- ✓ Good purity with simple purification.
- ✓ Good productivity, ~g/L.



-Template, Temperature, Time-

ver. 2 (4 mM GSH), 1 ng/μL DNA
 reaction at 16/23/30/37°C for 4/8/24 hours

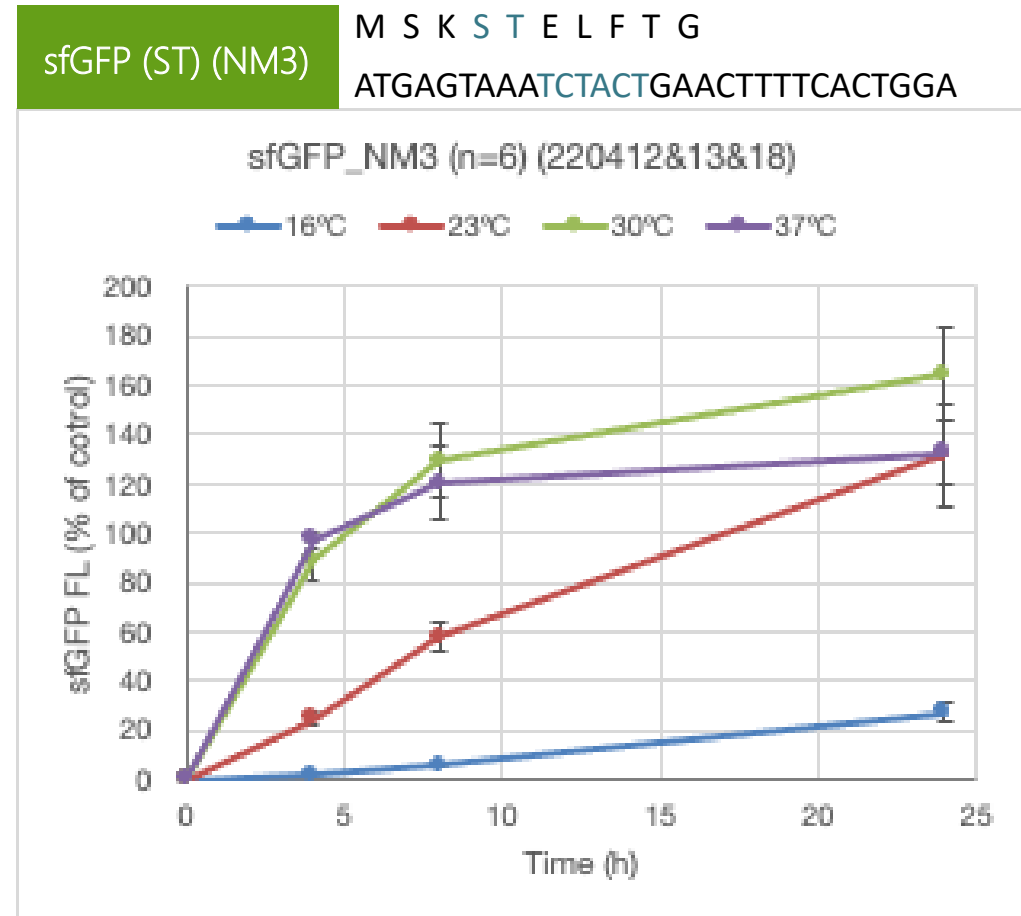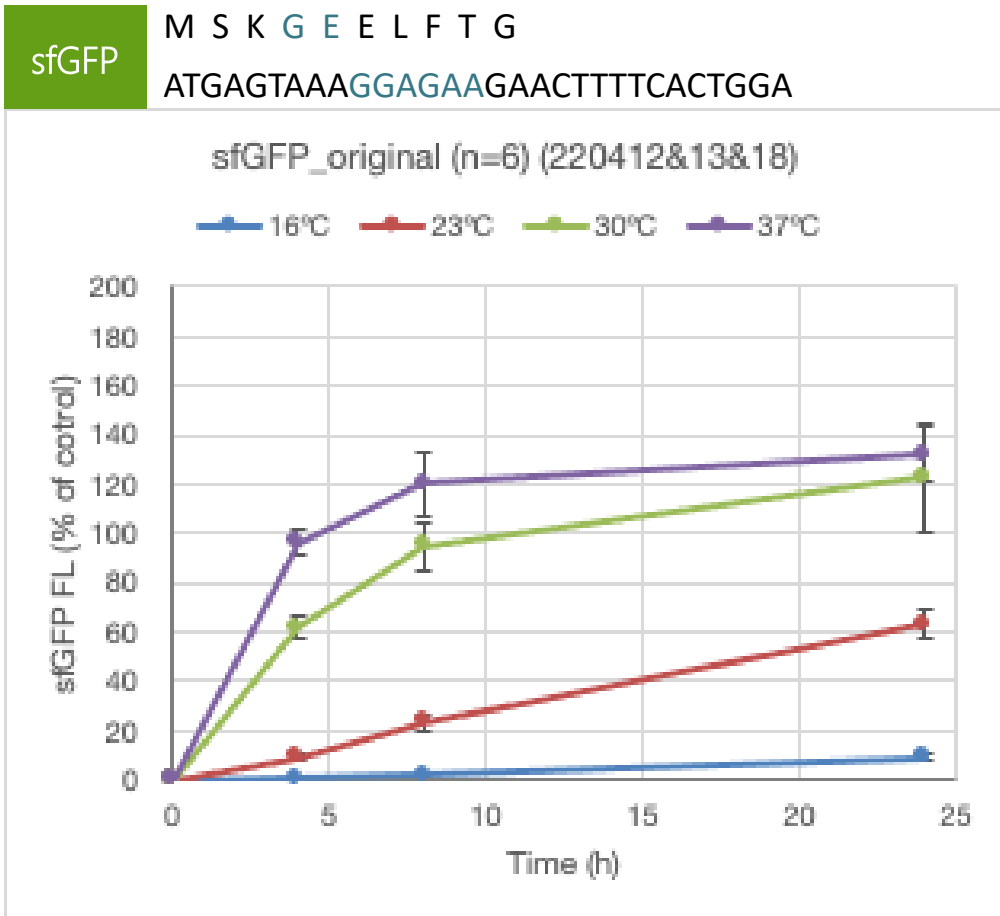




PUREfres[®]


-Improve process from Days to Hours-

PUREfres[®] Synthesis

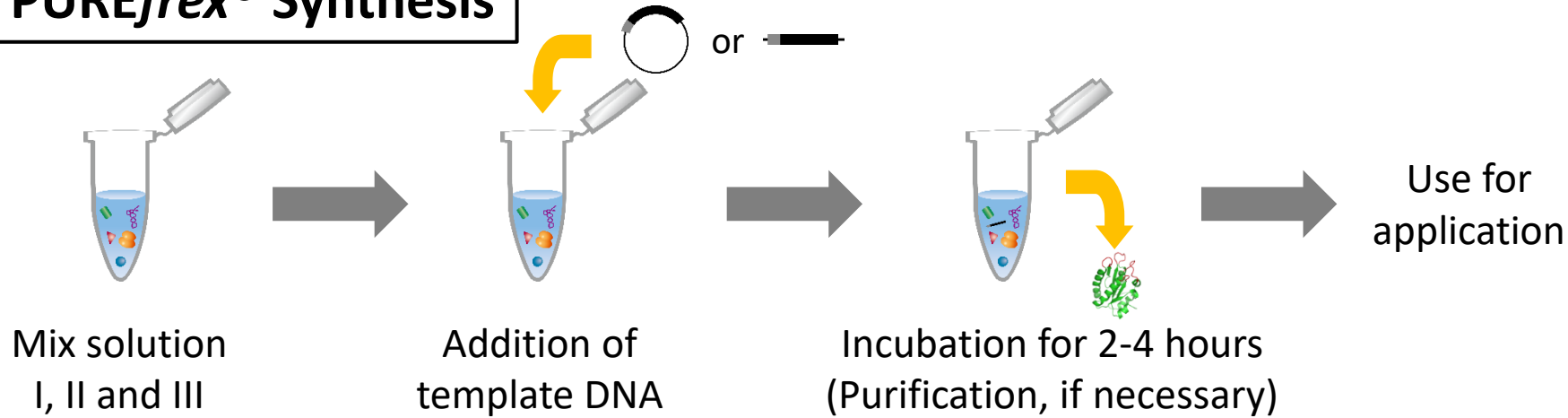

Total: 2-4 hours


E. coli Expression

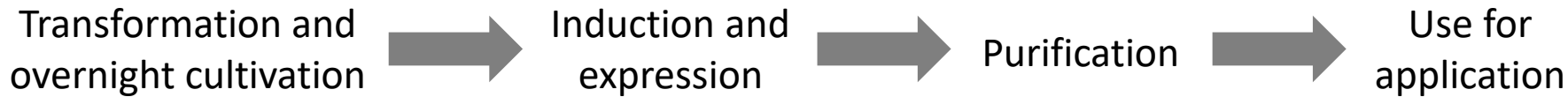


Total: 3-4 days

Design
Suitable for highspeed learning

Express → Check

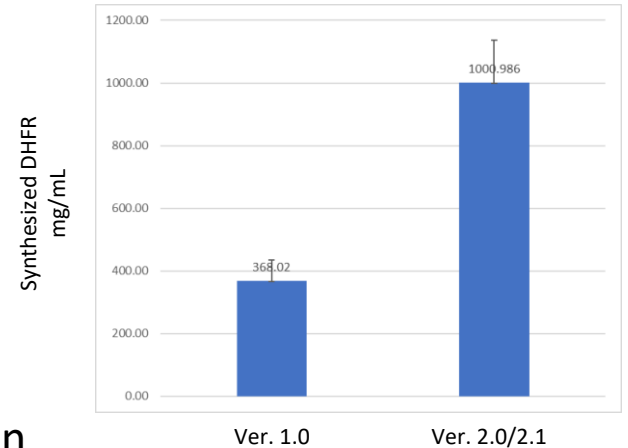



μL ~ *L scale*

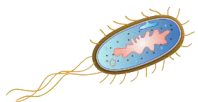




-Expression of scFv and Fab-


SDS-PAGE

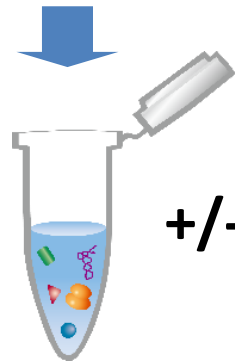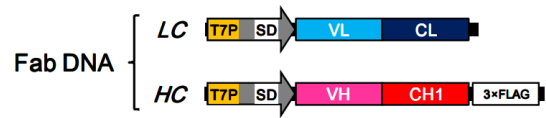

+/- **DsbC Set**
DnaK Mix
For correct folding

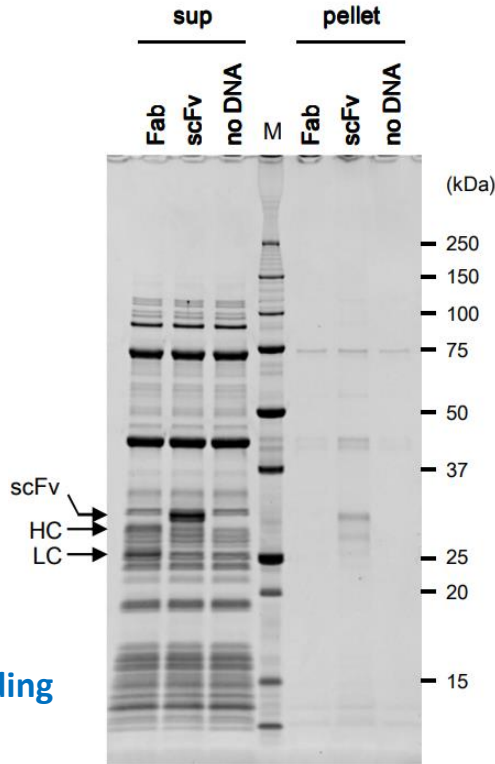

Activity

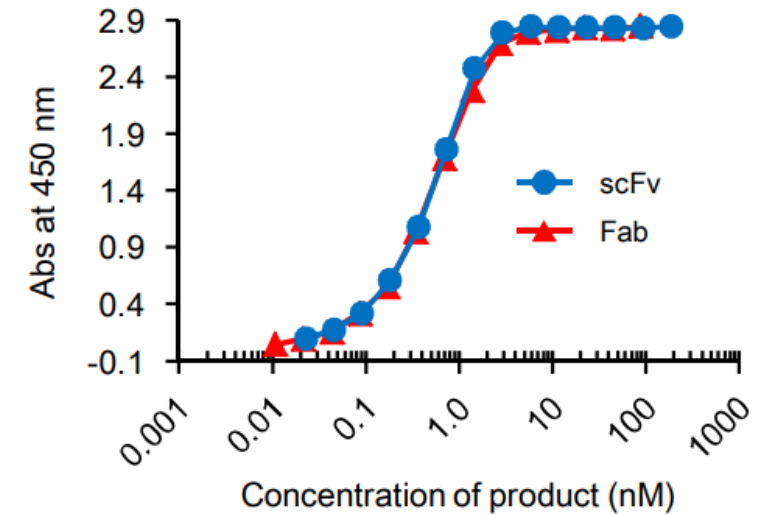


[Murakami et al. \(2019\) Sci. Rep. vol.9, p.671. \(Supplementary Information\)](#)




✓ **Active scFv and Fab (derived from Herceptin) can be synthesized.**



In vitro expression and Biacore analysis of Fab fragments

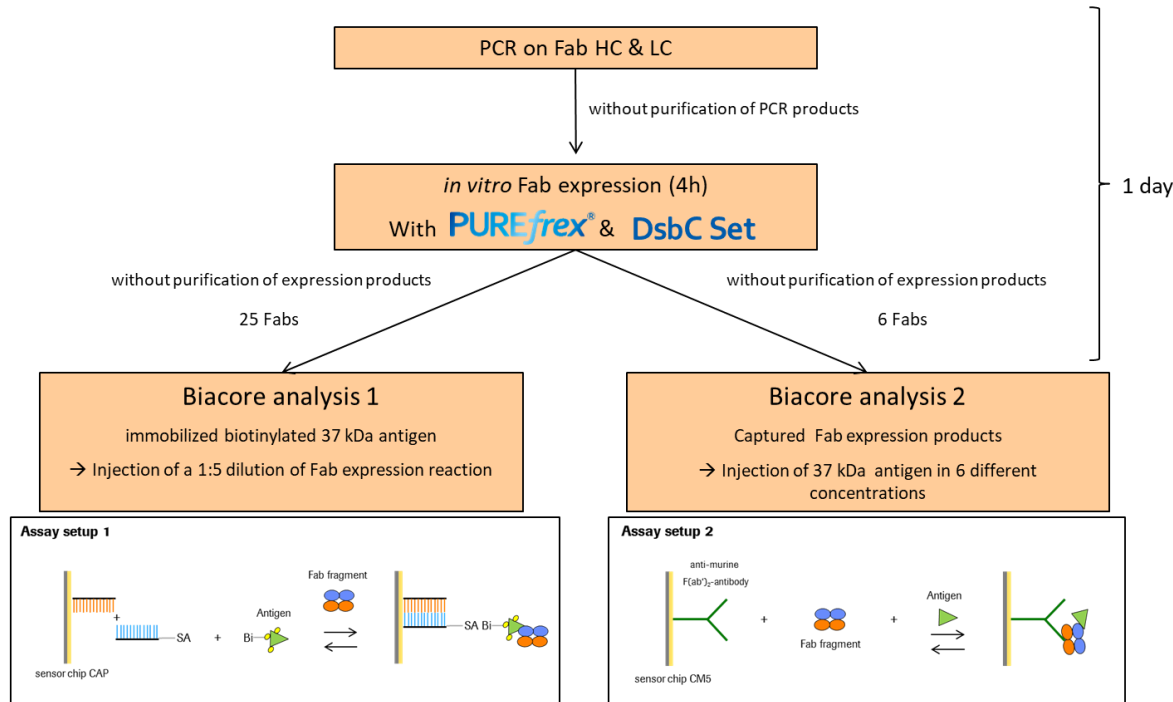

Kinetic analysis of 25 Fab binders

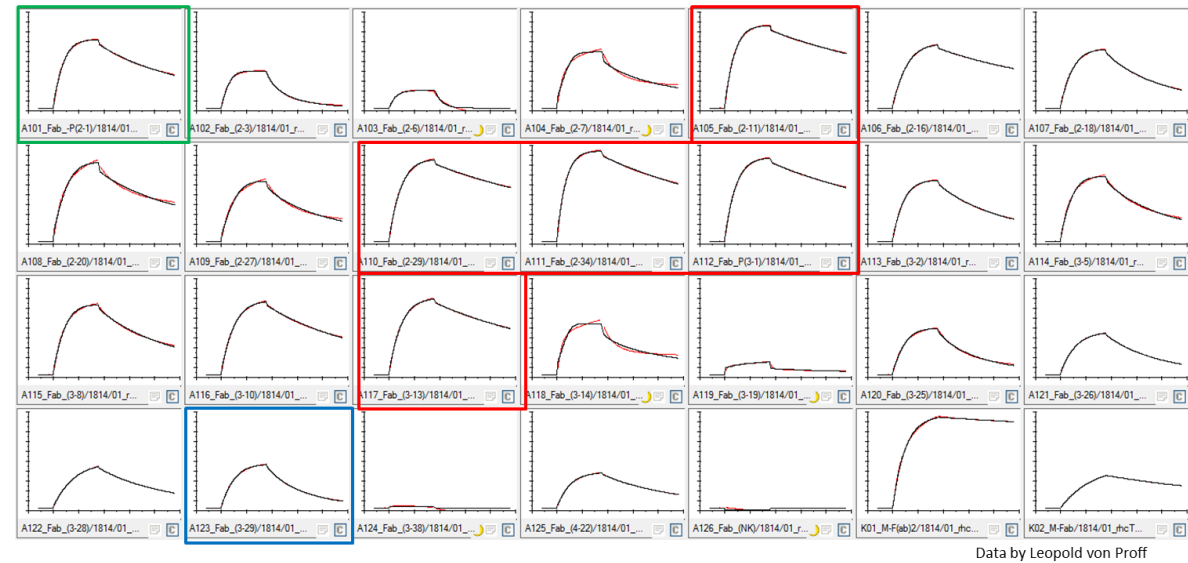


→ Selection of Fabs for further kinetic analysis

✓ **Active Fabs are expressed/screened in HT manner.**



iBody's Ecobody Technology

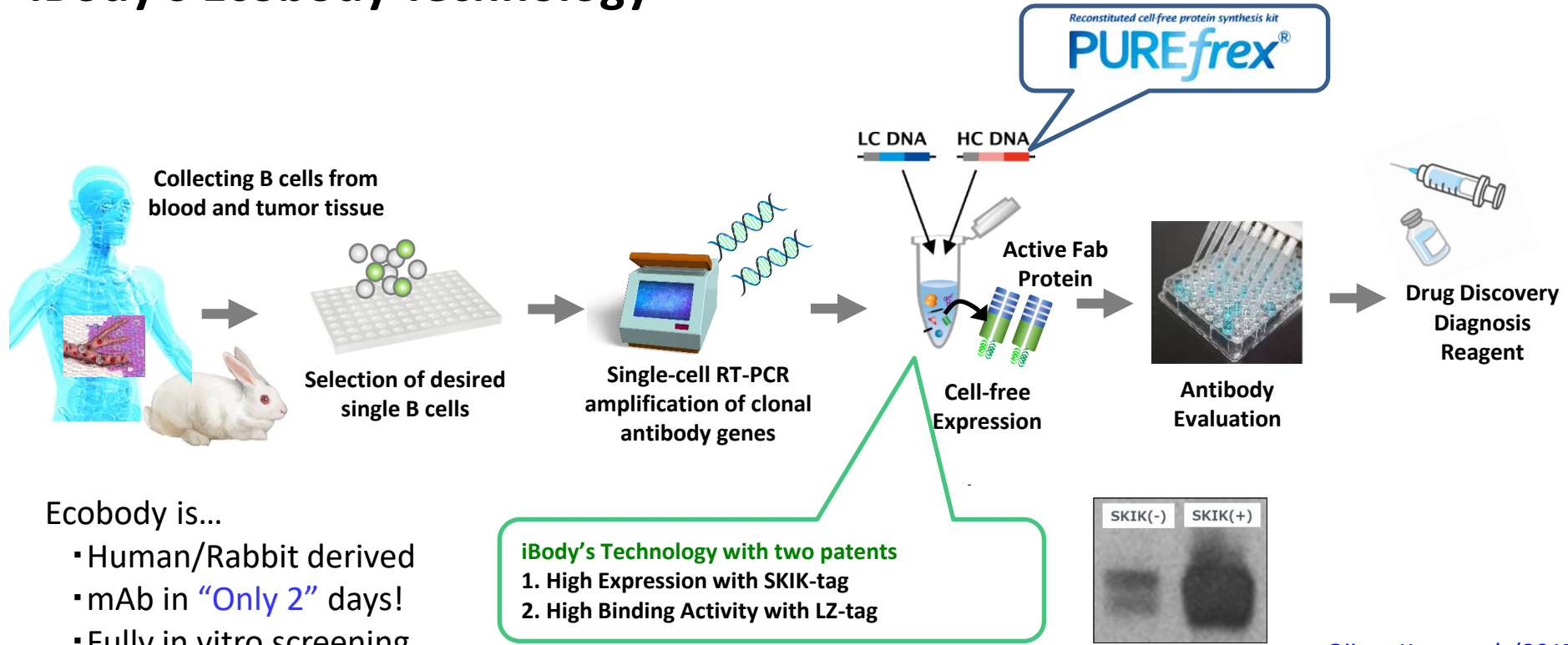

Ecobody is...

- Human/Rabbit derived
- mAb in "Only 2" days!
- Fully in vitro screening
- No culture

[Ojima-Kato et al. \(2017\) Sci. Rep., 7, 13979.](#)

<https://www.ibody.co.jp/en/>




✓ Active Fab is expressed/screened in HT manner.



SDNC system

(*In vitro* Selection by Differencing operation using NGS and Cell-Sorting with cDNA display)

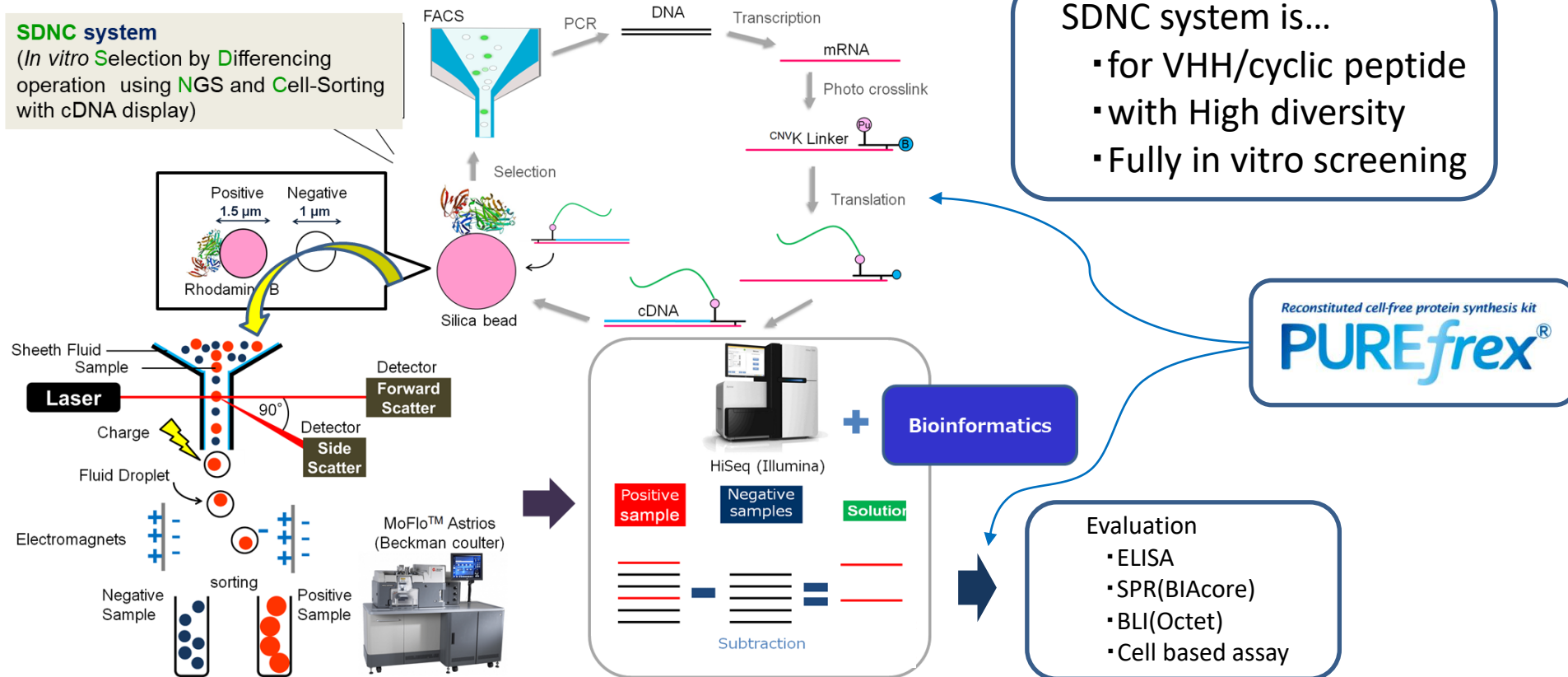


<https://www.epsilon-mol.co.jp/eng/>

✓ PUREfres is applied for cDNA display based screening.







-Application for BiTE-

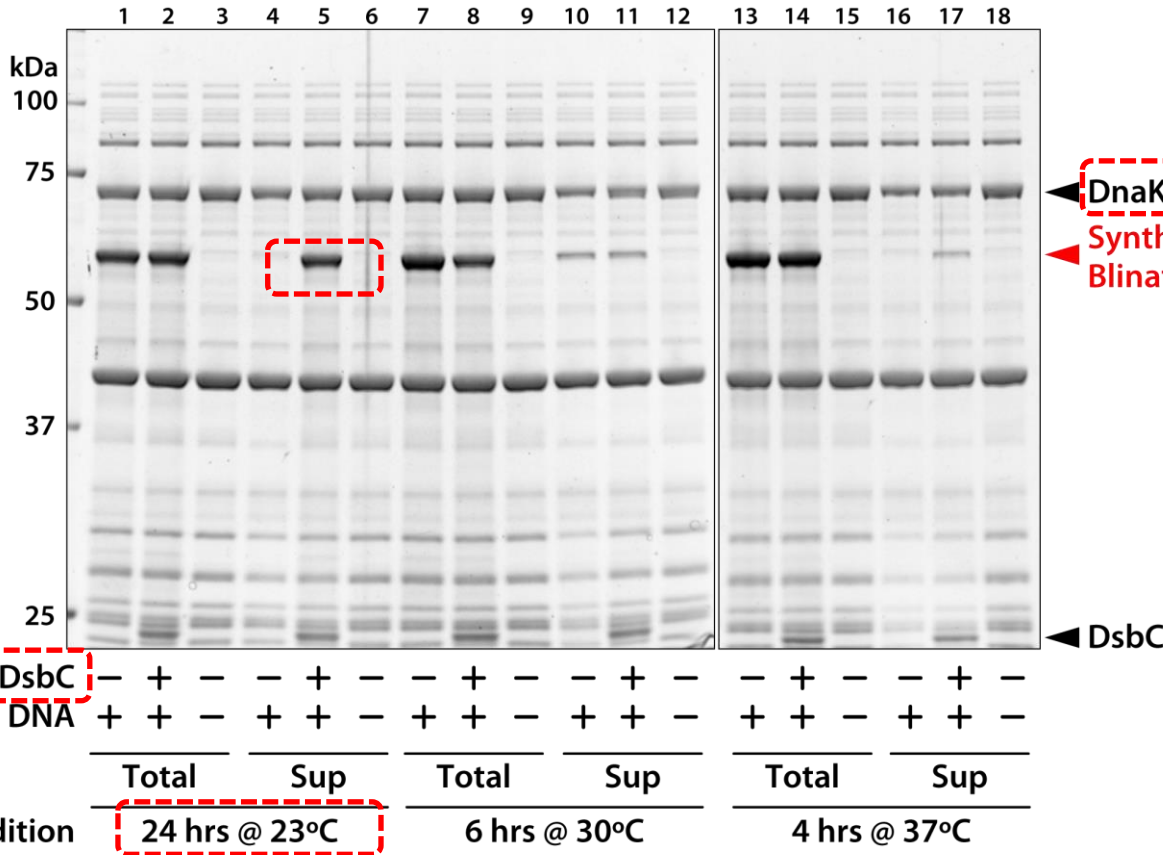


anti-CD3-scFv/CD19-scFv

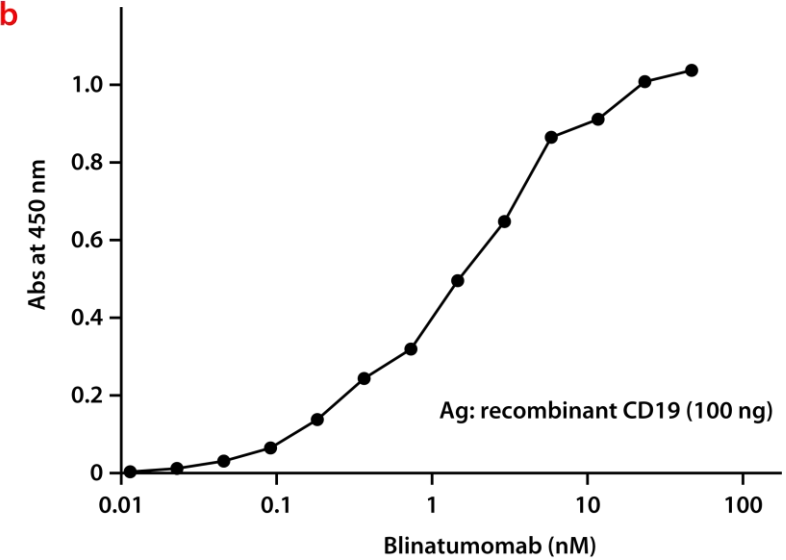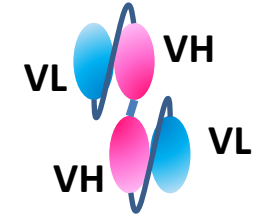



✓ 0.6mg/mL active BiTE was produced under optimized condition.



-Application for IgG-

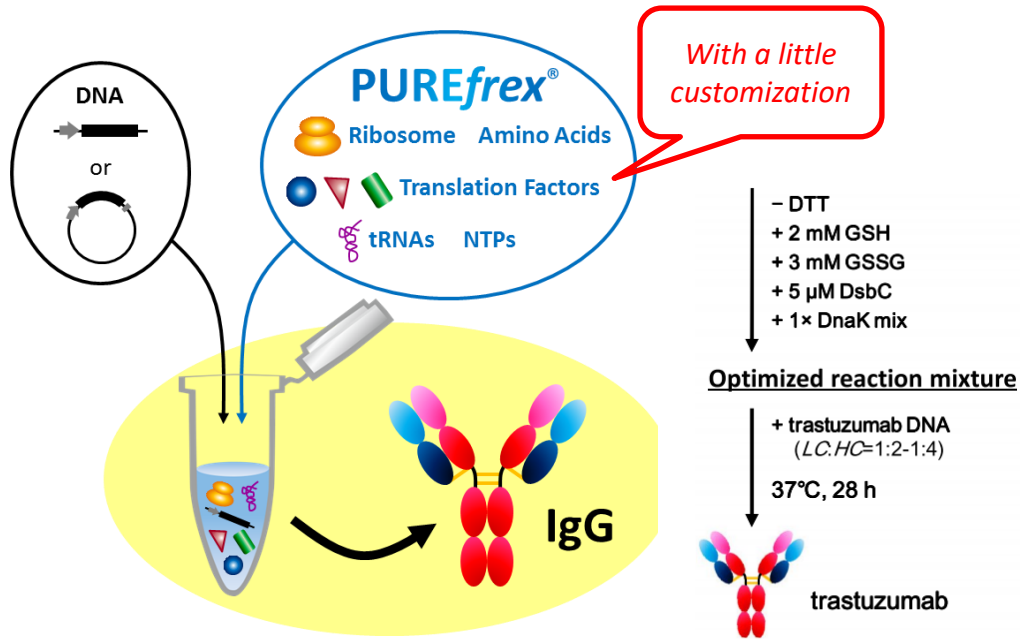

With a little customization

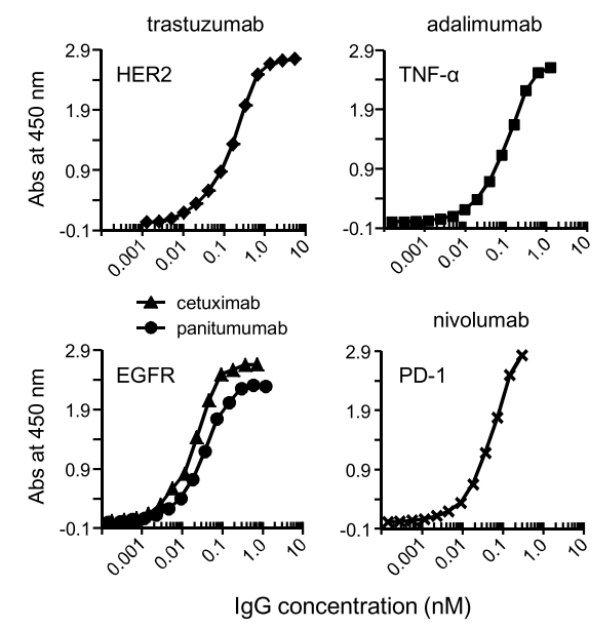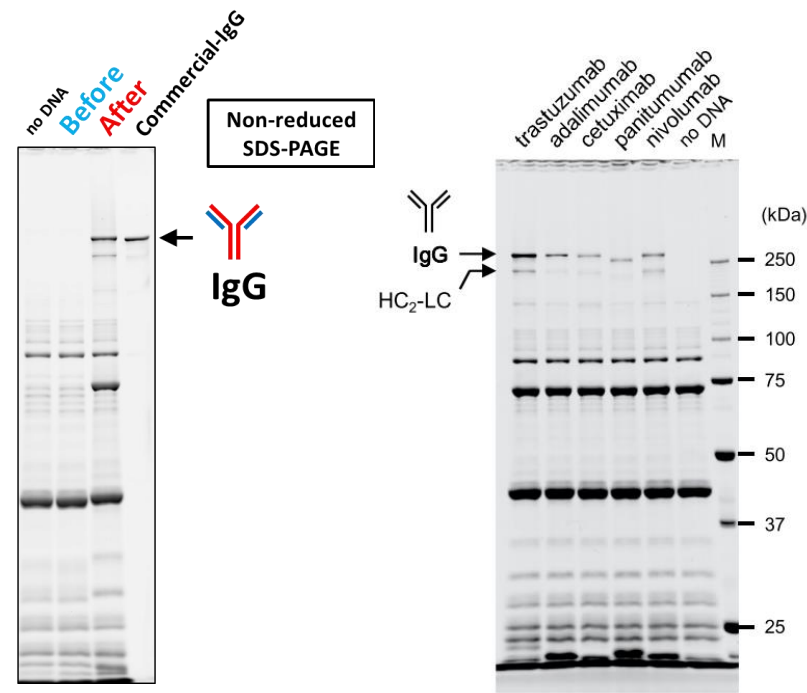

Only 2 days for IgG!!



[Murakami et al. \(2019\) Sci. Rep. vol.9, p.671.](#)

✓ Full size IgG can be synthesized.







-Application for complex molecule-

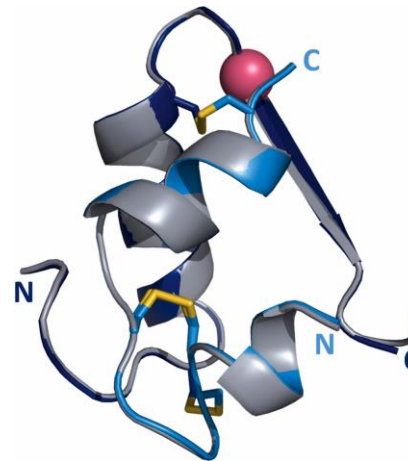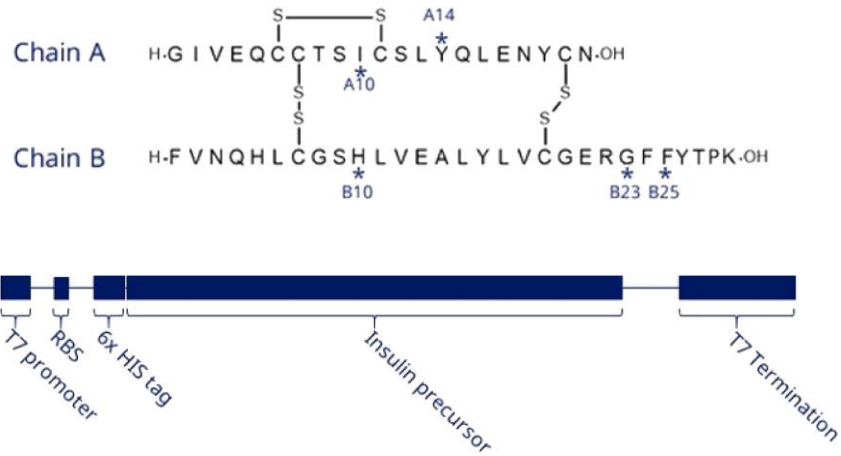

-	-	+	-	-	+	-	-	+	-	-	+	-	-	+	4 mM GSH/3 mM GSSG
-	+	+	-	+	+	-	+	+	-	+	+	-	+	+	0.375 mg/ml DsbC
+	+	-	+	+	-	+	+	-	+	+	-	+	+	-	2 mM DTT

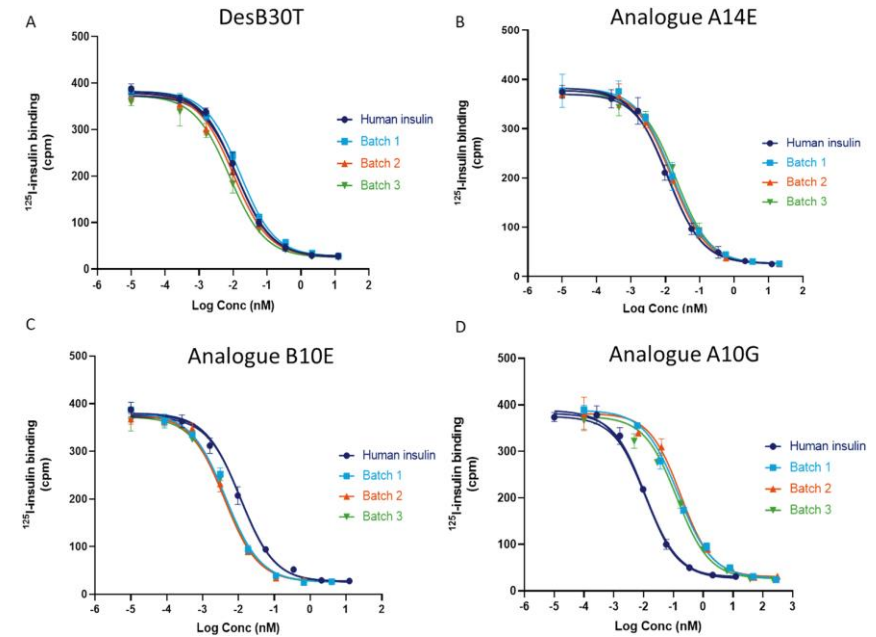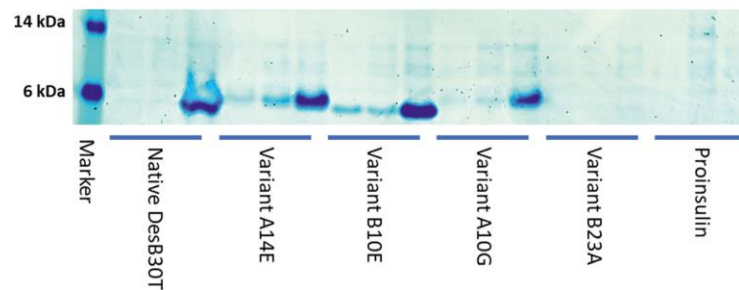


[Jensen et al. \(2021\) Protein Expr. Purif., 186, 105910.](#)



✓ **Functional insulin analogues can be synthesized.**




-Application for glycosylation; CFPS-IVG-

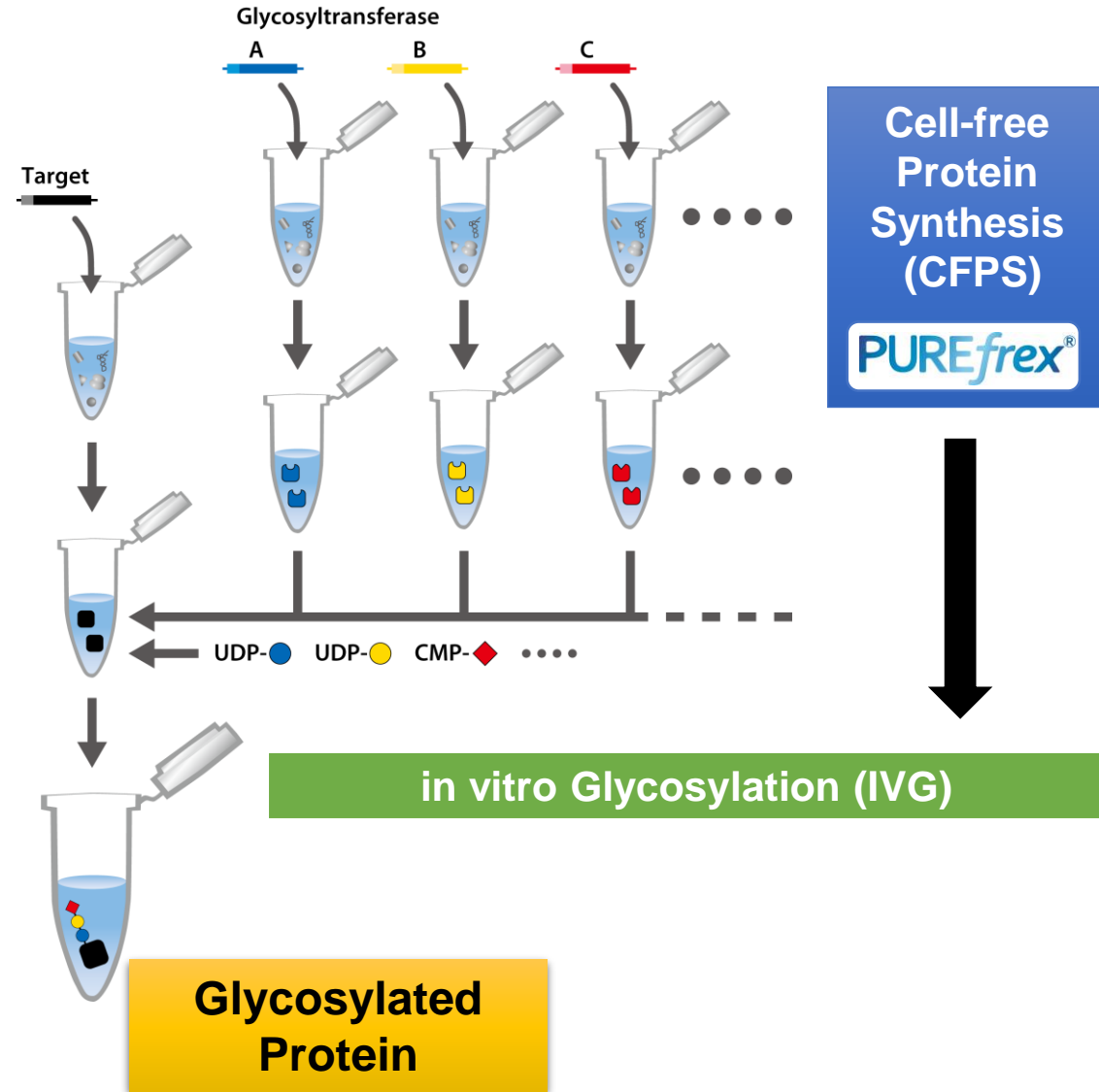

(kDa)

15

10

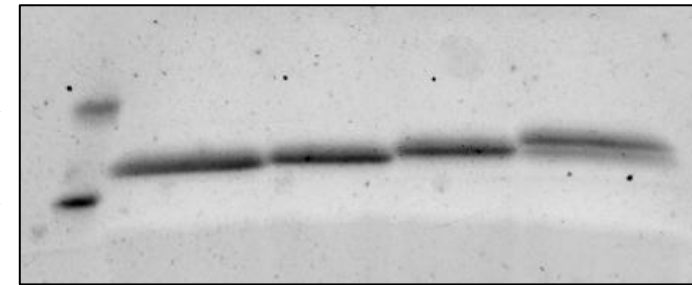





UDP-● / Enzyme A	-	+	+	+
UDP-● / Enzyme B	-	-	+	+
CMP-◆ / Enzyme C	-	-	-	+

**Applicable to glycosylation of target proteins.
Various reaction system can be designed.**





-Application for glycosylation; CFPS-IVG-

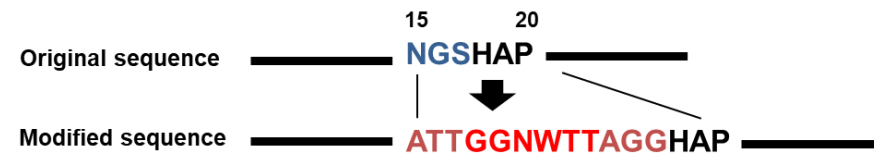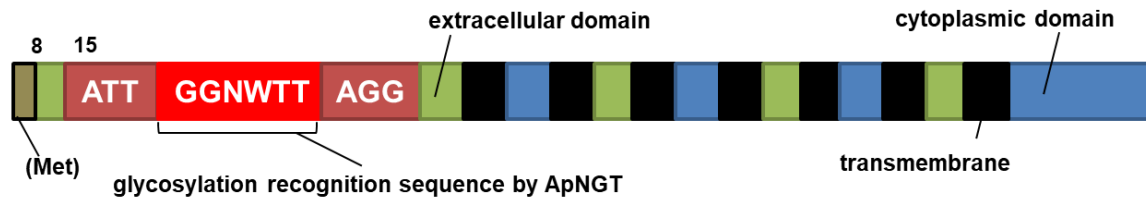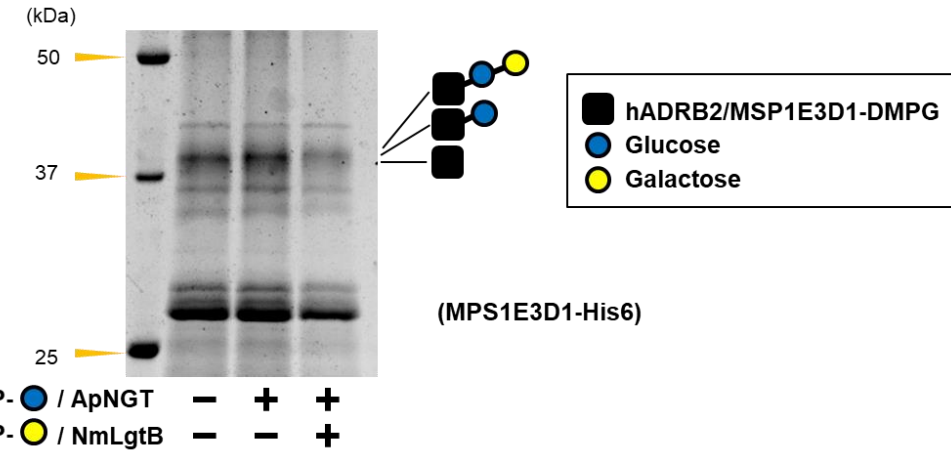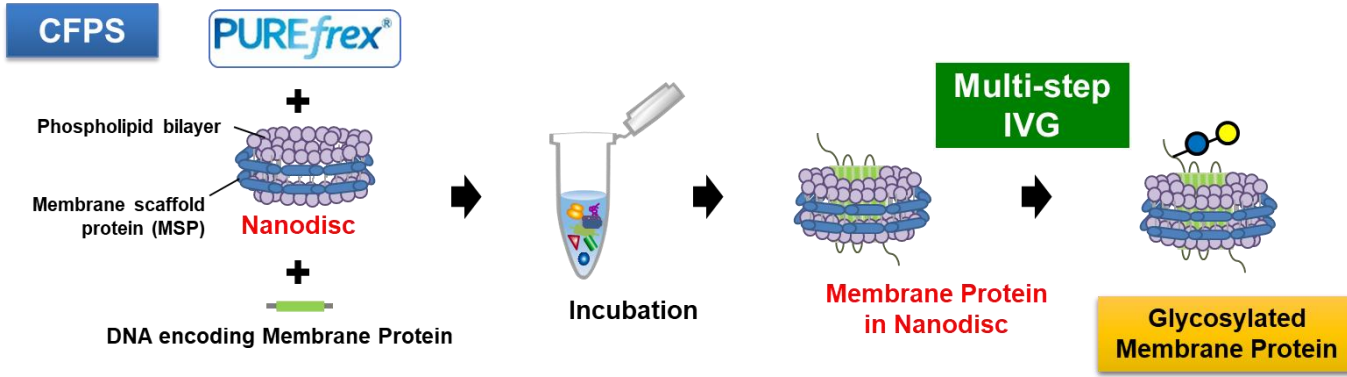

Solubilized membrane protein, hADRB2, are also glycosylated by multi-step IVG






-Application for acetylation-


Acetylation (by NatB)

target: α -synuclein (K6A)

MDVFMAGLSK... Hisx6

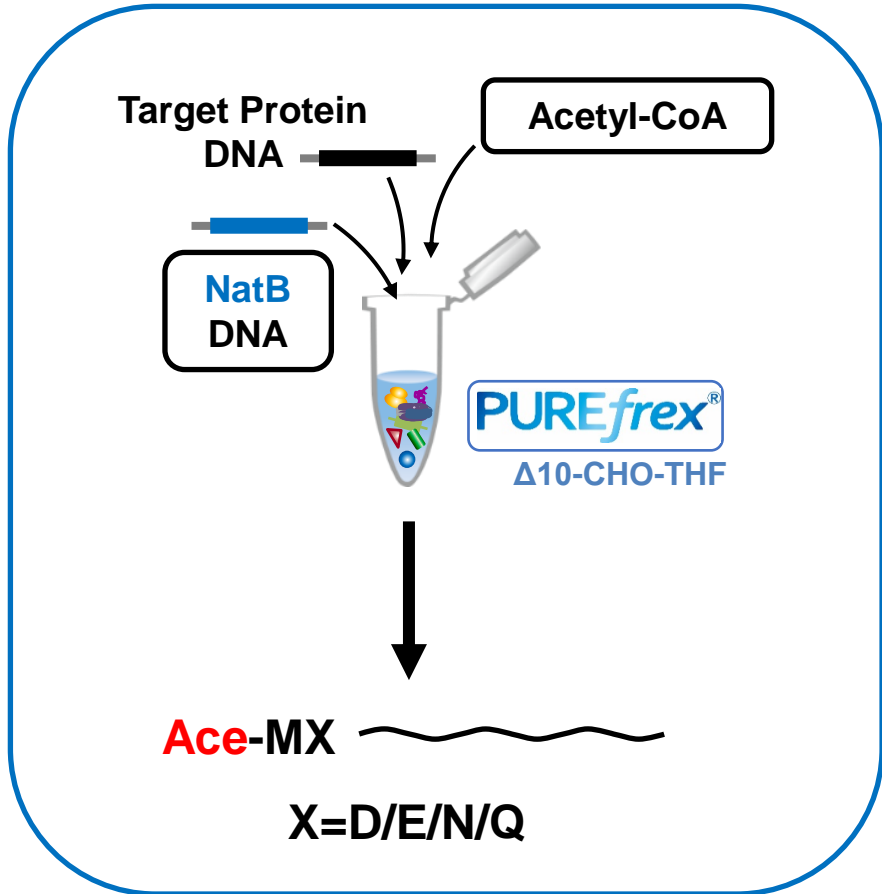

Chromatogram of MS1 intensity

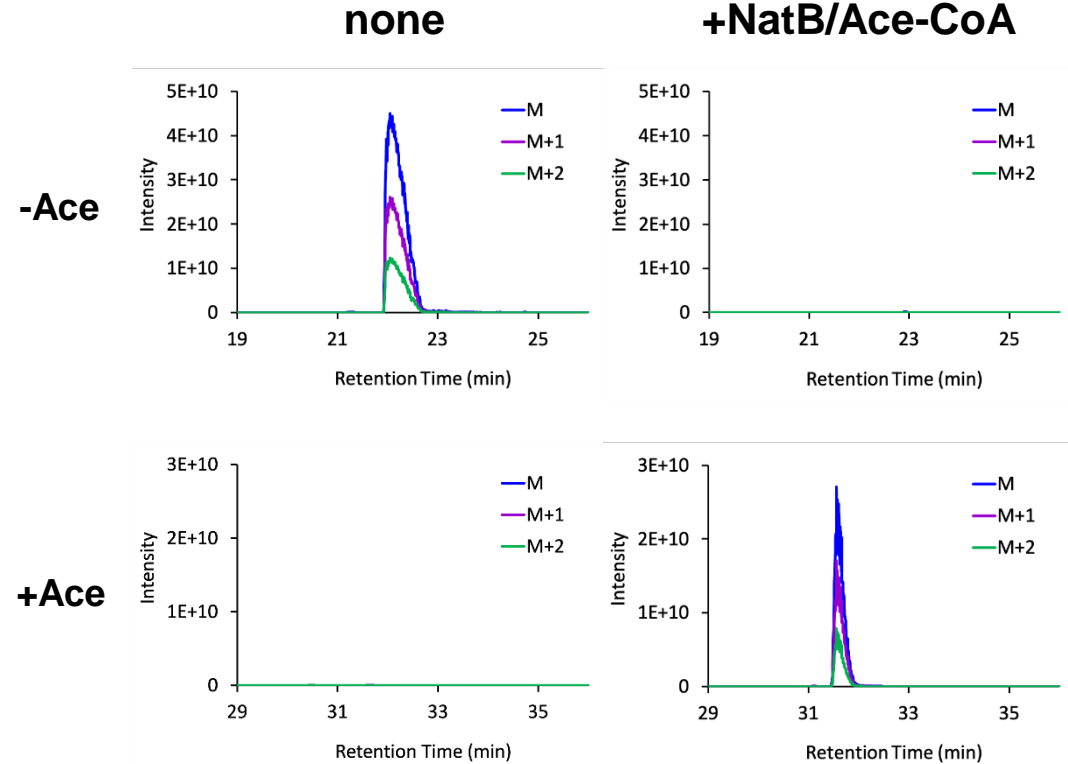

Perfect control of N-terminal acetylation of α -synuclein




Ribosome display system using PUREfres

◆ Advanced screening system for Biologics


- mAb (scFv / Fab)
- VHH
- Cyclic peptide

◆ High Selection Efficiency

- Completely molecular based system
- >10¹² diversity

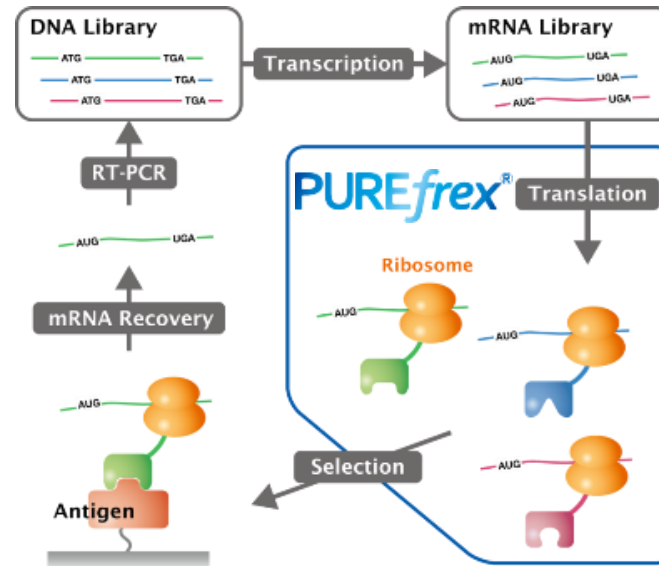

Licensed technology under JP4931135 etc.

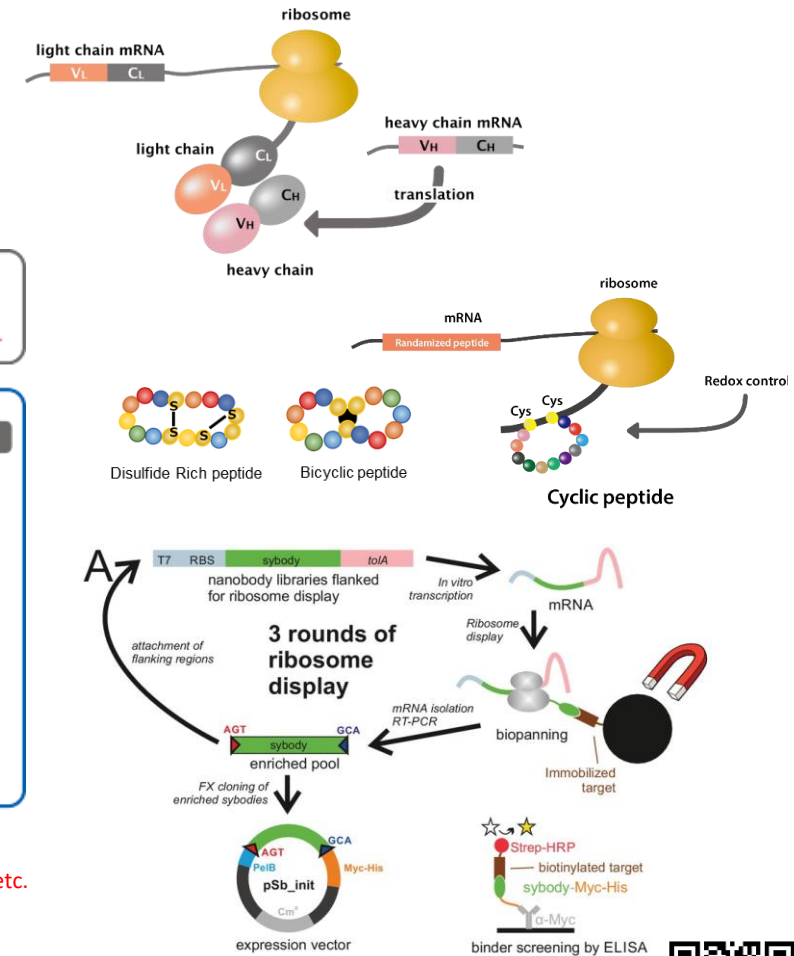


Zimmermann I. et al. (2018) eLife, 7, e34317.







-Broad applications, yet to come!-

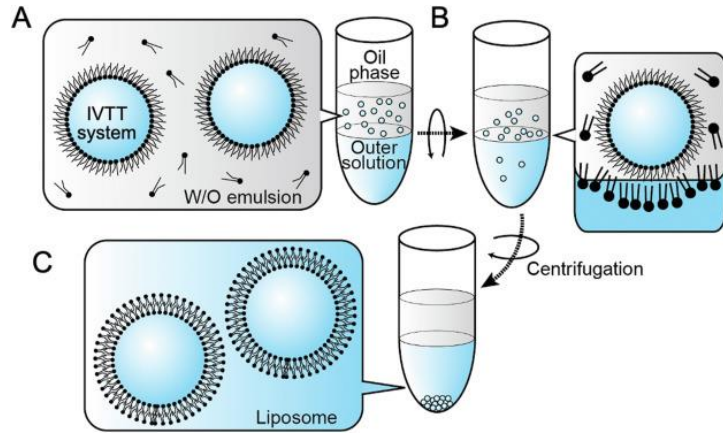


[Uyeda et al. \(2022\) J. Biosci & Bioeng. vol.133, p181.](#)

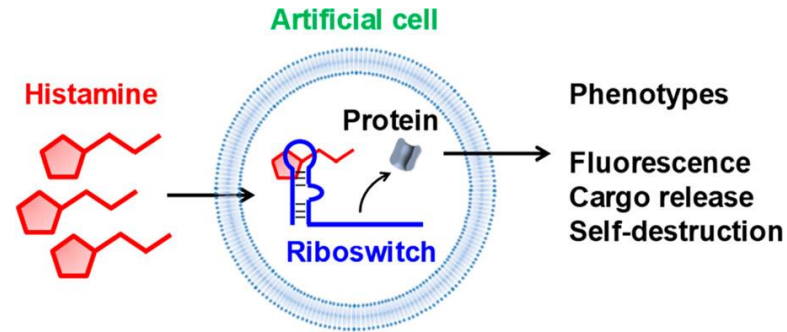


[Dwidar et al. \(2019\) J. Am. Chem. Soc. vol.141, p.11103.](#)

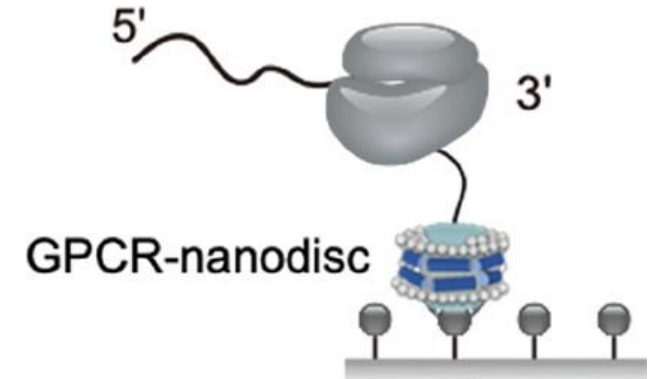


[Nakai et al. \(2022\) Anal. Chem. vol.94, p.3831.](#)

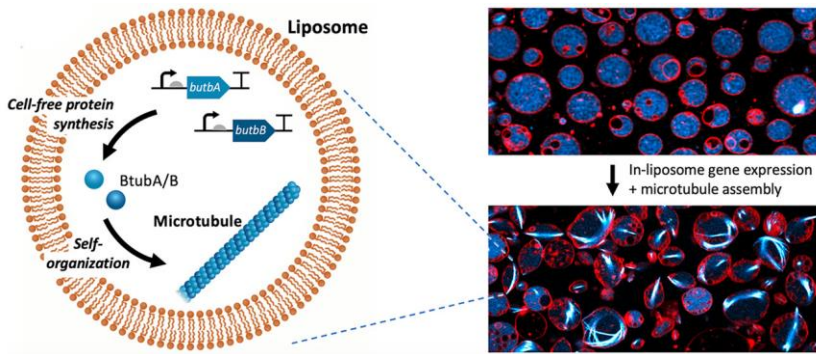

[Kattan et al. \(2021\) ACS. Synth. Biol. vol.10, p.2447.](#)

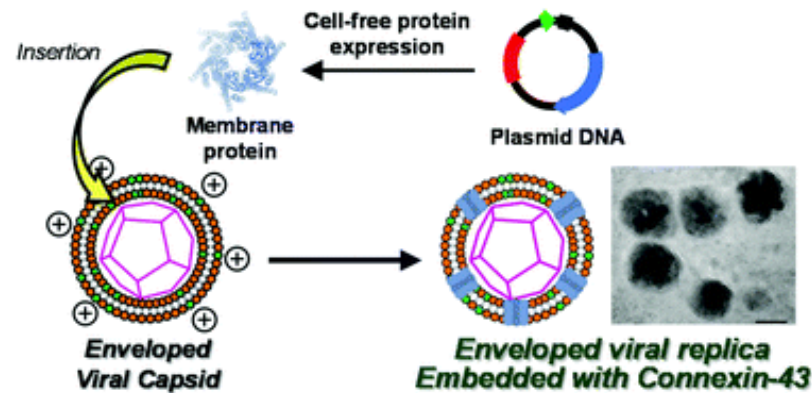

[Furukawa et al. \(2022\) RSC. Chem. Biol. vol.3, p.231.](#)

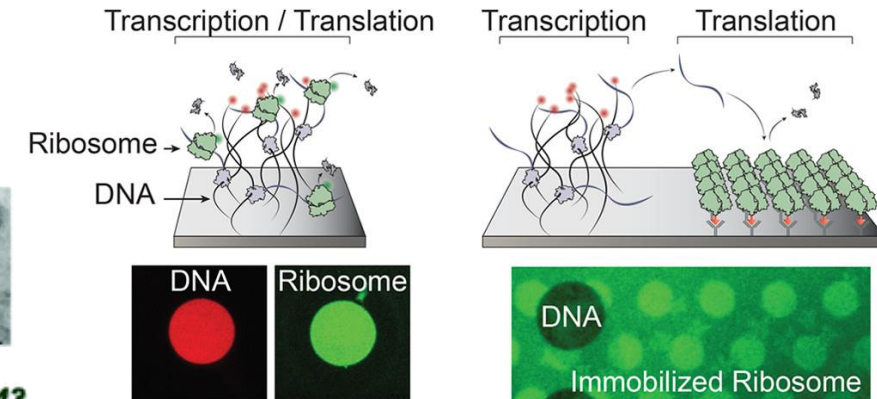

[Levy et al. \(2021\) ACS. Synth. Biol. vol.10, p.609.](#)

Contact information

Booth# 61



Reconstituted cell-free protein synthesis kit

PUREfrex[®]

For reagent use for expression and screening of biologics

<https://purefrex.genefrontier.com/?pegs2022-eu.html>



in vitro protein selection technology

PUREfrex[®]RD

*For screening service / collaboration / technology transfer
for generation of new biologics*

Takashi Ebihara, Ph.D., COO, GeneFrontier

E-mail: ebihara@genefrontier.com