

PUREfrex® 2.1

#PF213-2ML-EX For 2 mL Reaction
 #PF213-10ML-EX For 2 mL x 5 Reaction
 #PF213-50ML-EX For 2 mL x 25 Reaction

In vitro research use only
 Store at -80°C before opening

Introduction

About PUREfrex®

PUREfrex® is a reconstituted cell-free protein synthesis kit which GeneFrontier has developed based on the PURE system technology. The target protein can be synthesized by adding the template DNA (or mRNA) to the reaction mixture. The PURE system is a unique cell-free protein synthesis system, which has originally developed by Professor Takuya Ueda at the University of Tokyo, and consists of only purified factors necessary for transcription, translation and energy regeneration (Ref. 1). Therefore it enables to adjust the composition of the reaction mixture.

PUREfrex® has been raised in purity by improving the methods for preparing ribosomes, tRNAs and all proteins in the reaction mixture compared with the original PURE system (Ref. 2). As the result, the contaminating lipopolysaccharide from *E. coli* is reduced to less than 1 EU per 1 µL of reaction and other contaminants, such as RNase and β-galactosidase, are also reduced.

Because all of proteins in PUREfrex® have no tags, the synthesized protein can be purified and detected by any tags.

- (References) 1. Shimizu *et al.* (2001) *Nat. Biotechnol.*, vol. 19, p. 751
 2. Shimizu *et al.* (2005) *Methods*, vol. 36, p. 299

Kit components

	Volume	2ML	10ML	50ML
• Solution I *1	800 µL	×1	×5	×25
	Amino acids, NTPs, tRNAs and substrates of enzymes Store at -20°C			
• Solution II *2	100 µL	×1	×5	×25
	Proteins in 30% glycerol buffer Store at -20°C or -80°C *3			
• Solution III *2	200 µL	×1	×5	×25
	Ribosomes (20 µM) Store at -80°C *3			
• Cysteine *1	160 µL	×1	×5	×25
	Cysteine (10 mM) Store at -20°C			
• DTT *1	160 µL	×1	×5	×25
	Dithiothreitol (40 mM) Store at -20°C			
• GSH *1	160 µL	×1	×5	×25
	Reduced glutathione (80 mM) Store at -20°C			
• DHFR DNA *2, 4	10 µL	×1	×1	×1
	PCR product (20 ng/µL) containing a gene encoding <i>E. coli</i> DHFR Store at -20°C			

Store at -80°C before opening

*1) Reducing agents are excluded from Solution I in PUREfrex® 2.1. You can freely adjust the amount of reducing agents (Cysteine, DTT and GSH, etc.), **but be sure to add Cysteine, which is a material of protein.**

The standard working concentration of Cysteine, DTT and GSH are 0.5, 2 and 4 mM, respectively. Other reducing agents are also available for the reaction, e.g. 2-mercaptoethanol and TCEP.

We recommend optimizing the concentration of reducing agent to synthesize a protein with higher activity because the suitable concentration is depends on the synthesized protein.

*2) Solution II, Solution III and DHFR DNA are the same as PUREfrex® 2.0 (#PF201).

*3) For storage at -80°C, the remaining solution should be frozen rapidly in liquid nitrogen or a freezing mixture of dry ice and ethanol. Please divide into aliquots, if necessary, and avoid refreeze and thaw as much as possible.


*4) As a positive control for the protein synthesis reaction, 1 µL of DHFR DNA should be added to 20 µL of reaction. Please visit our web site to get the nucleic acid sequence of DHFR DNA.

How to aliquot

For instructions on how to aliquot each component, please refer to our website or the "PUREfrex® 2.1 Instruction Manual."

Template DNA

1. Construction of the template DNA

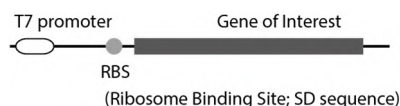

The template DNA for the protein synthesis by PUREfrex® should contain T7 promoter and ribosome binding site (SD sequence) at the upstream of the gene encoding the target protein.

At the downstream of stop codon, more than 10 nucleotides are required.

All three stop codons are available.


Both circular and linear DNAs are available as a template DNA.

For the circular DNA, e.g. plasmid DNA, T7 terminator is required at the downstream of stop codon.

For the linear DNA, e.g. PCR product or DNA fragment digested by restriction enzymes, T7 terminator is not necessarily required.

2. Preparation of the template DNA

An example of the method for preparing the template DNA by two-step PCR is shown in below.

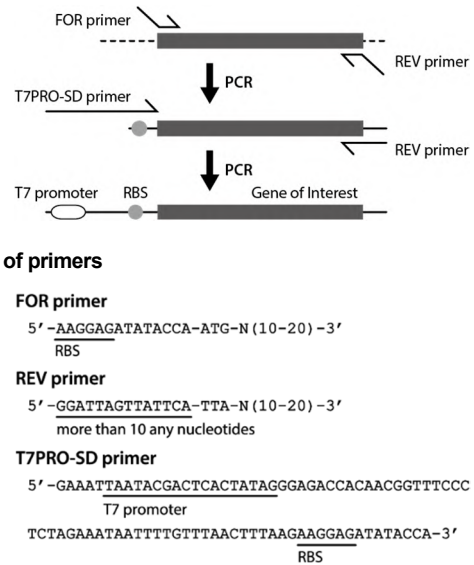

3. Sequence of primers

FOR primer

5' -AAGGAGATATACCA-ATG-N (10-20) -3'
RBS

REV primer

5' -GGATTAGTTATTC-ATA-N (10-20) -3'
more than 10 any nucleotides

T7PRO-SD primer

5' -GAAATTAATACGACTCACTATAGGGAGACCAACGGTTTCCC
T7 promoter
TCTAGAAATAATTTTGTTTAACTTTAAGAAGGAGATATACCA-3'
RBS

Protocol

The protein synthesis reaction using PUREfrex® 2.1 can be performed in any volume. For example, 20 µL of reaction mixture containing 0.5 mM Cysteine and 4 mM GSH is assembled as below.

1. Thaw completely Solution I, Cysteine and GSH by incubation at room temperature or 37°C for 5 minutes for completely dissolving, and then leave at room temperature.
2. Thaw Solution II and Solution III on ice.
3. Mix each solution by vortex and centrifuge briefly to collect solution at the bottom.
4. Assemble the reaction mixture in a tube as follows.
(Add the template DNA to 0.5-3 ng/µL per 1 kbp)

Water	7-X	µL
Solution I	8	µL *5
Cysteine	1	µL
GSH	1	µL
Solution II	1	µL
Solution III	2	µL
Template DNA	X	µL
Total	20	µL

5. Incubate the tube at 37°C for 4-6 hours.
6. Analyze the synthesized product.

*5) Please note that the volume of Solution I in PUREfrex® 2.1 (#PF213) is different from PUREfrex® 2.0 (#PF201).

Note

PUREfrex® is developed for *in vitro* research use only. PUREfrex® should not be used for the therapy, diagnostic or administration to animals including human and should not be used as food or cosmetics etc.

To avoid the contamination of nuclease, nuclease-free-treated water, reagents and materials should be used. We also recommend wearing gloves and mask.

PUREfrex® is a registered trademark of GeneFrontier Corporation. For any commercial use of PUREfrex®, please contact GeneFrontier in advance.

e-mail : purefrex@genefrontier.com

Distributor



COSMO BIO USA

【Outside Japan】
 2792 Loker Ave West, Suite 101, Carlsbad, CA 92010, USA
 email: info@cosmobiousa.com
 Phone/FAX: (+1) 760-431-4600
 URL: www.cosmobiousa.com



COSMO BIO CO., LTD.

【JAPAN】
 TOYO EKIMAE BLDG. 2-20, TOYO 2-CHOME,
 KOTO-KU, TOKYO 135-0016, JAPAN
 Phone: +81-3-5632-9610 FAX: +81-3-5632-9619
 URL: https://www.cosmobio.co.jp/

Manufacturer



GeneFrontier
 www.genefrontier.com

GeneFrontier Corporation
 SHARP Kashiwa Building, 4F, 273-1 Kashiwa,
 Kashiwa-shi, Chiba 277-0005 Japan